


## LabGenius Mini BlockHeater Operating Manual

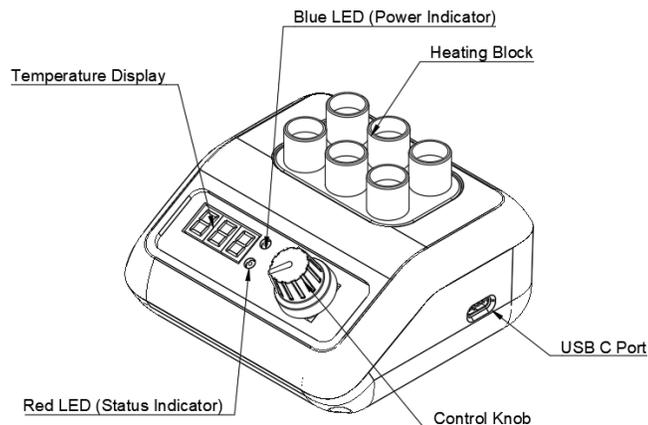


### Set up:

Power the BlockHeater with a 5V USB C power adapter. The LED display will turn on promptly and indicate the current temperature value in °C.

To change the temperature set up, press the control knob one time till the red LED is blinking. Rotate the knob to adjust the set up temperature. Press the control knob again till the red LED is solid and wait till temperature reach its set up temperature.

To change the heating block, simply lift up the current well plate and insert the desire well plate.

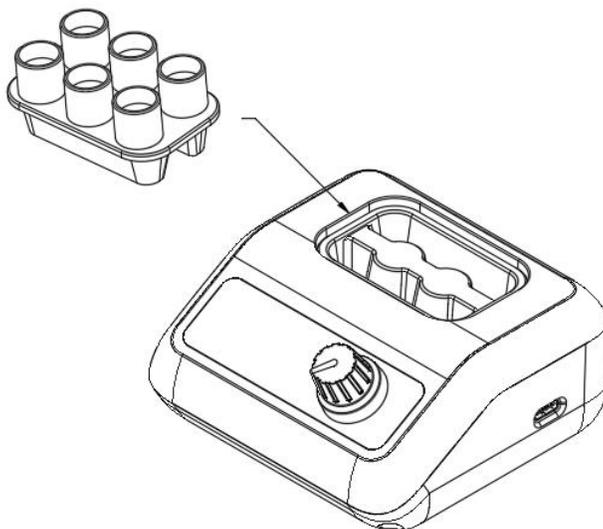

**Note:** Please do not touch the heating block while temperature is hot. Touching the heating block while hot can cause severe burn and injury to the operator.