Instruction Manual

Corning® Scholar™ 170 Hot Plate and Scholar™ 171 Stirrer

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Plaque Chauffante Corning Scholar 170 et Agitateur Scholar 171

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Corning® Scholar™ 170 Hot Plate and Scholar 171 Stirrer
(100V, 120V, 230V)

Contents

1.0	International Symbols	2
2.0	Product Description	2
3.0	Operating Conditions	2
4.0	Warnings	2
5.0	Heating Operation and Controls	3
6.0	Stirring Operation and Controls	3
7.0	Inspection Procedure	3
8.0	Preventive Maintenance	4
9.0	Technical Specifications	4
10.0	Technical Support/Service	4
11.0	Warranty Registration	4
12.0	Warranty	5
13.0	Product Disposal	6

1.0 International Symbols

Hot surface indicator, surfaces on this unit may be hot to the touch.

This unit may cause the emission of toxic gasses due to heating of certain materials.

Earth ground

2.0 Product Description

Model	Top Style	Top Size (in.)	Dimensions (in.)
Corning® Scholar™ 170 Hot Plate	Enameled	5 x 5	4 x 5 x 5
Corning Scholar 171 Stirrer	Plastic	5 x 5	3.2 x 5 x 5

3.0 Operating Conditions

The Corning Scholar is designed to be safe when used under the following conditions:

- ▶ Indoor use
- ▶ Altitude up to 2,000 meters
- ▶ Temperature of 0°C to 40°C
- ▶ Maximum relative humidity of 80% for temperatures up to 31°C decreasing linearly to 50% relative humidity at 40°C.
- ▶ Position at least 12" from walls, 48" from ceiling, and 12" between hot plates if using multiple units.
- ▶ Main supply voltage fluctuations not to exceed $\pm 10\%$ of the nominal voltage.
- ▶ Installation Category II
- ▶ Pollution Degree 2

4.0 Warnings

- ▶ **DO NOT** immerse unit for cleaning.
- ▶ **DO NOT** remove or modify ground power plug. Use only properly grounded outlets to avoid shock hazard.
- ▶ **DO NOT** use insulating materials, such as a sand bath, on top of this product when heating.
- ▶ **USE CAUTION** when heating volatile materials, top surface and element can reach the "Flash Point" temperature of many chemicals.

- ▶ **NOT EXPLOSION PROOF OR SPARK PROOF.**

- ▶ **USE** appropriate hand and eye protection when handling hazardous chemicals.
- ▶ **DO NOT** modify the unit electrically or mechanically without Corning's written approval as personal injury or product damage may occur.

5.0 Heating Operation and Controls

The Corning® Scholar™ 170 Hot Plate will not heat when the dial setting is turned to the "OFF" position. To operate the unit turn the dial until the dial indicator lines up with the desired dial setting. The following table details approximate temperatures for various dial settings.

Dial Setting	Approx. Temperature
Low	60°C
2	100°C
3	140°C
4	160°C
5	200°C
6	230°C
High	360°C

NOTE: Temperatures shown are approximate "no load" values. Variables affecting actual temperature may include environment conditions, dial setting position, and top loading.

6.0 Stirring Operation and Controls

The Corning Scholar 171 Stirrer will not stir when the dial setting is turned to the "OFF" position. To operate the unit, turn the dial until the dial indicator lines up with the desired dial setting. The stir range is 100 to 1,600 rpm.

7.0 Inspection Procedure

- ▶ **Look for etching or chipping on top plate.** If any of these conditions are present, discontinue use and contact Corning for additional instructions.
- ▶ **Look for damaged power cord.** If the cord is damaged, do not modify the grounded power plug to avoid shock hazard or any other unit component. Service should always be done by qualified technicians using Corning instructions. Contact Corning for additional instructions.

8.0 Preventive Maintenance

- ▶ **Disconnect power** by unplugging unit before performing any maintenance or service.
- ▶ **Always use a grounded power outlet.**
- ▶ **Keep top surface clean and dry.** Unplug unit and use non-abrasive cleaner. Contact Corning if there is any doubt about the compatibility of decontamination or cleaning agents.
- ▶ **Clean spills promptly.** If liquid or solid material gets inside the product, immediately disconnect power to the product, use non-abrasive cleaner, and discontinue use. Contact Corning for additional instructions regarding interior spills.

9.0 Technical Specifications

Corning Cat. No.	Model	Power	Temp. Range	Stir Range (rpm)
6795-170	PC-170 Hot Plate	120V/50-60Hz/ 250W/2.1A	60°C - 360°C	N/A
6796-170	PC-170 Hot Plate	230V/50-60Hz/ 250W/1.1A	60°C - 360°C	N/A
6797-170	PC-170 Hot Plate	100V/50-60Hz/ 250W/2.5A	60°C - 360°C	N/A
6795-171	PC-171 Stirrer	120V/50-60Hz/ 30W/0.25A	N/A	100 - 1,600

10.0 Technical Support/Service

If you have a question or a service inquiry about the Corning® Scholar™ 170 Hot Plate or Scholar 171 Stirrer, contact Corning Customer Service immediately at 1.800.492.1110 (United States and Canada), 1.978.442.2200 (outside the United States), or Contact your local Corning sales office.

Before returning any unit for service, a Return Authorization (RA) number must first be obtained. Equipment sent in without our prior authorization will be returned at the customer's expense. When returning a unit to Corning for service, it should be sent in the original packaging. If this is not possible, be sure that the unit is sufficiently packed. Any damage resulting from improper packaging is the responsibility of the customer. A written explanation should accompany the unit along with the RA number.

11.0 Warranty Registration

Your purchase of this product must be registered with Corning in order for your warranty to be in effect. Please register your warranty online at www.corning.com/lifesciences/warranty.

12.0 Warranty

Corning Incorporated (Corning) warrants that this product will be free from defects in material and workmanship for a period of (1) year from date of purchase. CORNING DISCLAIMS ALL OTHER WARRANTIES WHETHER EXPRESSED OR IMPLIED, INCLUDING ANY IMPLIED WARRANTIES OF MERCHANTABILITY OR OF FITNESS FOR A PARTICULAR PURPOSE. Corning's sole obligation shall be to repair or replace, at its option, any product or part thereof that proves defective in material or workmanship within the warranty period, provided the purchaser notifies Corning of any such defect. Corning is not liable for any incidental or consequential damages, commercial loss or any other damages from the use of this product.

This warranty is valid only if the product is used for its intended purpose and within the guidelines specified in the supplied instruction manual. This warranty does not cover damage caused by accident, neglect, misuse, improper service, natural forces, or other causes not arising from defects in original material or workmanship. This warranty does not cover motor brushes, fuses, light bulbs, batteries, or damage to paint or finish. Claims for transit damage should be filed with the transportation carrier.

In the event this product fails within the specified period of time because of a defect in material or workmanship, contact Corning Customer Service at the following numbers: USA: 1.800.492.1110; Canada: 1.978.442.2200. For other regions of the world, visit www.corning.com/lifesciences or see the included instruction manual for a list of worldwide support offices.

Corning's Customer Service team will help arrange local service where available, coordinate a return authorization number, or provide a return authorization number and shipping instructions. Products received without proper authorization will be returned. All items returned for service should be sent postage prepaid in the original packaging or other suitable carton, padded to avoid damage. Corning will not be responsible for damage incurred by improper packaging. Corning may elect for onsite service for larger equipment.

Some states do not allow limitation on the length of implied warranties or the exclusion or limitation of incidental or consequential damages. This warranty gives you specific legal rights. You may have other rights which vary from state to state.

No individual may accept for, or on behalf of Corning, any other obligation of liability, or extend the period of this warranty.

For your reference, make a note of the model and serial number, date of purchase, and supplier here.

Model No. _____ Serial No. _____

Date Purchased _____

Supplier _____

13.0 Product Disposal

According to Directive 2012/19/EU of the European Parliament and Council of 4th July 2012 on waste and electronic equipment (WEEE) as amended, Corning® Scholar™ 170 Hot Plate or Scholar 171 Stirrer is marked with the crossed-out wheeled bin and must not be disposed of with domestic waste.

Consequently, the buyer shall follow the instructions for reuse and recycling of waste electronic and electrical equipment (WEEE) provided with the products and available at the following link: www.corning.com/weee.

Register your product warranty online at www.corning.com/lifesciences/warranty.

For more specific information on claims, visit the Certificates page at www.corning.com/lifesciences.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

For additional product or technical information, visit www.corning.com/lifesciences or call 800.492.1110. Outside the United States, call +1.978.442.2200 or contact your local Corning sales office.

Corning Incorporated Life Sciences

836 North St.
Building 300, Suite 3401
Tewksbury, MA 01876
t 800.492.1110
t 978.442.2200
f 978.442.2476

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For a listing of trademarks, visit www.corning.com/clstrademarks.
All other trademarks are the property of their respective owners.

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